

# GUARDIAN CLEAN

## Machine Warewashing Tablets

---

**Guardian Clean Tablets** are suitable for all warewasher-safe glassware. The product formula especially protects glasses against corrosion and clouding. Combined with outstanding cleaning performance, Guardian Clean dissolves heavy deposits such as protein and starch build-up. The tablet application is designed for easy and economic use.

**Guardian Clean** has been shown under Laboratory conditions to assist in achieving clean and sanitary conditions of your ware when used at temperatures of 55°C and higher as required by your Safety Programme.



## WAREWASHER TABLETS

**Guardian Clean Tablets** are designed for manual and machine cleaning of laboratory glassware

**Guardian Clean Tablet** is a non-caustic, non-chlorine, non-phosphate formula with biodegradable surfactants to remove even the toughest soils in one pass

**Guardian Clean** has been shown to assist with reducing commonly occurring pathogens *Escherichia coli* and *Listeria monocytogenes*. Under laboratory conditions >log 5 reductions were observed when **Guardian Clean** was used at 55°C for a 20 minute contact time. Use concentration = 1 tab per 8.5 litres of water, typifying domestic-style warewasher conditions.

NB: Test result available on request

### DIRECTIONS FOR USE

- ▲ Check product label for detailed use information
- ▲ Take one tablet of **Guardian Clean Tablet** and open the plastic foil
- ▲ Place the tablet in the product reservoir and start normal program
- ▲ For large machines, dose directly into the wash tank



### SAFETY INFORMATION

Do not mix with any other products. For further information and safety instructions see product label and Material Safety Data Sheet.

#### For professional use only

This detergent is ideally suited to under bench machines. Select low temperature washing cycles, wash full loads and do not exceed the recommended dosage. This will minimise both energy and water consumption without compromising sanitation result.

Recommended storage temperature between 0°C and 40°C in original container only.

Product Code	Product Description	Pack Size
19046400	Guardian Clean Machine Warewashing Tablets	4kg / 200 Tablets

PLEASE SPEAK TO YOUR HEALTHCARE MANAGER OR CALL **1800 022 002** FOR FURTHER INFORMATION